For Research Use Only. Not for use in diagnostic procedures.

V5-tagged Protein PURIFICATION KIT

Ver.2

(MoAb. clone OZA3)

CODE No.3317

*PURIFICATION to Maintain Protein Activity
from eukaryote cell lysate and culture supernatant*

MBL MEDICAL & BIOLOGICAL LABORATORIES CO., LTD.

URL <http://ruo.mbl.co.jp>

e-mail support@mbi.co.jp, TEL 052-238-1904

Product Description

The ability to isolate and study a purified protein lies at the heart of modern biochemistry. Researchers in many fields require highly purified, active proteins for studies involving signaling pathways, enzymology, receptor binding, DNA binding, post-transcriptional modifications, and much more. Thus, choosing a method of purification is an important aspect in maintaining protein structure and function.

Recombinant tagged protein purification methods and kits are now widely recognized. The V5 epitope tag, GKPIP NLLGLDST, is derived from the P/V proteins of paramyxovirus SV5. This V5-tag is also commonly used in such as mammalian and insect cell expression vectors.

MBL's V5-tagged Protein PURIFICATION KIT Ver.2 is designed for the isolation of V5-tagged protein from cell culture supernatants containing serum and cell lysate under neutral pH condition. Severe conditions such as acidic or alkaline elution denature protein structure. However, a neutral pH elution can preserve protein activity and native conformation. MBL has developed the Anti-V5-tag Beads to purify V5-tagged proteins quickly and efficiently. As the Beads can be used at neutral pH, the purified proteins can maintain the activity and conformation. The elution of V5-tagged proteins from the Beads is achieved by the addition of the V5 peptide. As the V5 peptide competes with V5-tagged proteins on the Beads, the purified proteins do not lose the protein activity. The simple procedures of this kit have been optimized by using a Spin Column resulting in high efficiency.

Kit Components

Components are sufficient for conducting 20 times purifications of V5-tagged protein.

- | | |
|-----------------------------|--|
| 1. Anti-V5-tag Beads | 25% slurry: 100 μ L beads in 400 μ L total volume in PBS with 0.1% ProClin 150 as preservative |
| 2. Elution Peptide | V5 peptide, 2 mg in 1 mL PBS after reconstitution |
| 3. Spin Columns Sets | 20 columns with pre-inserted bottom plugs and top caps |
| 4. Wash Concentrate | 10x concentrate, 6 mL |

Storage

Store for up to 1 year from date of receipt at 2-8°C. Do not freeze.

Product Capacity

The purification capacity of the Anti-V5-tag Beads varies depending upon the V5-tagged protein. For examples, 5 μ L of Anti-V5-tag Beads (20 μ L slurry) bound 8.2 μ g of a V5-tagged protein (32 kDa) and eluted 6.3 μ g of purified protein.

Materials Required but not Provided

1. Microcentrifuge capable of 15,000 x g
2. Sampling tube (1.5 mL)
3. End-over-end rotator
4. PBS
5. Lysis buffer

Suitable Lysis buffer varies with cell type.

Note: see Additional Information

Homemade Lysis buffer

20-50 mM	Tris-HCl, pH 7.5 or HEPES-KOH, pH 7.5
50-250 mM	NaCl
5 mM	EDTA
1%	NP-40 or Triton X-100

if necessary add Protease Inhibitor Cocktail
(e.g. SIGMA code P8340, PIERCE code 78415).

Protocols

The following protocols are for the isolation of V5-tagged proteins produced in a 100-mm cell culture dish. The expression level of the V5-tagged protein may vary. If necessary, adjust the volume of Anti-V5-tag Beads and Elution Peptide Solution proportionally.

Material Preparation

1. Wash Solution

Dilute Wash Concentrate with 9 times its volume of distilled water.


(e.g. Dilute 0.1 mL of Wash Concentrate with 0.9 mL of distilled water.)

For each Spin Column, prepare 1 mL of Wash Solution.

2. Elution Peptide Solution

Reconstitute the Elution Peptide with 1 mL of distilled water. If you want to store the reconstituted Elution Peptide, prepare appropriate aliquots (e.g. 45 μ L x 20 tubes) and store at -20°C. Repeated freezing and thawing is not recommended.

Procedure Summary (Purification from mammalian cultured cell lysate)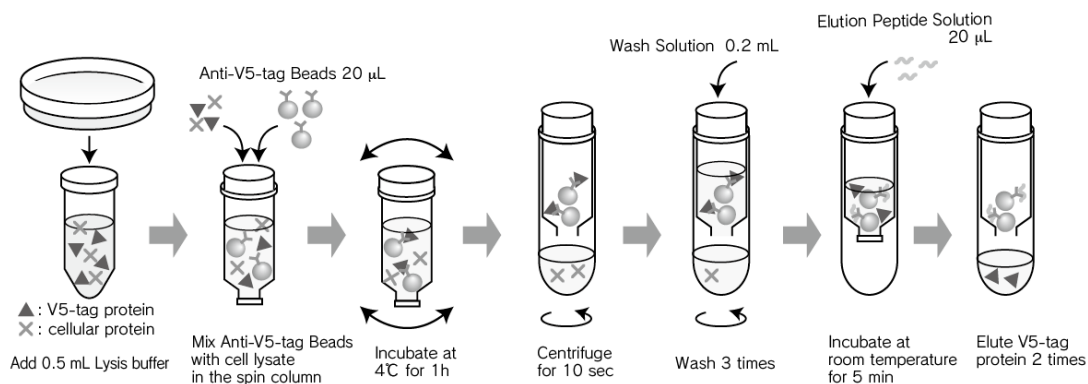

A. Purification from mammalian cultured cell lysate

(V5-tagged protein is not secreted from the cells)

(Lysis of Mammalian or Insect Cells)

1. Detach the cells from the culture dish if necessary, and collect the cell suspension into the centrifuge tube.
2. Centrifuge the cell suspension at 400 x g for 5 minutes to pellet the cells. Carefully remove and discard the supernatant.
3. Wash cells by resuspending the cell pellet in ice-cold PBS.
4. Centrifuge the cell suspension at 400 x g for 5 minutes to pellet the cells. Carefully remove and discard the supernatant.
5. Add 0.5 mL of Lysis buffer to the cell pellet and vortex.
6. Sonicate the sample for 15 seconds.
7. Incubate the sample for 15 minutes on ice.
8. Remove cell debris by centrifugation at 15,000 x g for 5 minutes at 4°C.

(Purification of V5-tagged Protein)

9. Transfer the 0.5 mL of cell lysate (supernatant from step 8) to the Spin Column.
10. Resuspend the Anti-V5-tag Beads by tapping and inverting the vial several times immediately before dispensing. Don't vortex.
11. Dispense 20 μ L of Anti-V5-tag Beads suspension (5 μ L Beads) into the Spin Column. Screw on the cap.
12. Incubate with gentle end-over-end mixing for 1 hour at 4°C. If the Spin Column does not fit your end-over-end rotator, put it in a suitable tube (e.g. 15 mL centrifuge tube) that fits your end-over-end rotator.
13. Loosen the top cap on the column. Remove the bottom plug. Don't discard the bottom plug. Place the Spin Column in a sampling tube. Centrifuge for 10 seconds. Discard the flow-through (or save for future analysis).
14. Take off the top cap. Keep the bottom plug off. Place the Spin Column in a sampling tube.
15. Add 0.2 mL of Wash Solution to each column. It is not necessary to stir the Spin Column. Centrifuge for 10 seconds. Discard the flow-through. Repeat this step two additional times.
16. Place the Spin Column in a new sampling tube.
17. For the first elution, screw the bottom plug on tightly. Add 20 μ L Elution Peptide Solution to the Anti-V5-tag Beads, then screw the top cap on tightly. Tap the tube gently several times. Incubate for 5 minutes at room temperature. Remove the top cap and bottom plug. Centrifuge for 10 seconds.

18. For the second elution, it is not necessary to place the bottom plugs and top cap on the Spin Column. Add 20 μ L Elution Peptide Solution to the Anti-V5-tag Beads, then tap the tube gently several times. Incubate for 1 minute at room temperature. Centrifuge for 10 seconds.

The two eluates (step 17 and 18) may be pooled in one sampling tube.

B. Purification from culture supernatant

(V5-tagged protein is secreted into the culture supernatant)

1. Collect the culture supernatant from the cell culture dish into a 15 mL centrifuge tube.
2. Centrifuge at 400 x g for 5 minutes to remove cell debris.
3. Transfer the supernatant to a new 15 mL centrifuge tube.
4. Resuspend the Anti-V5-tag Beads by tapping and inverting the vial several times immediately before dispensing. Don't vortex.
5. Dispense 20 μ L Anti-V5-tag Beads suspension (5 μ L beads) into the 15 mL centrifuge tube with culture supernatant. Screw on the cap.
6. Incubate with gentle end-over-end mixing for 1 hour at 4°C.
7. Centrifuge at 400 x g for 5 minutes. Discard the supernatant (or save for future analysis), but leave 100-400 μ L supernatant above the Anti-V5-tag Beads.
8. Resuspend the Anti-V5-tag Beads in the 100-400 μ L supernatant by pipetting up and down several times.
9. Transfer the resuspended Anti-V5-tag Beads in 100-400 μ L supernatant to the Spin Column.
10. Keep the top cap off. Remove the bottom plug. Don't discard the bottom plug. Place the Spin Column in a sampling tube. Centrifuge for 10 seconds. Discard the flow-through (or save for future analysis). Keep the bottom plug off. Place the Spin Column in a sampling tube.
11. Add 0.2 mL of Wash Solution to each column. It is not necessary to stir the Spin Column. Centrifuge for 10 seconds. Discard the flow-through. Repeat this step two additional times.
12. Place the Spin Column in a new sampling tube.
13. For the first elution, screw the bottom plug on tightly. Add 20 μ L Elution Peptide Solution to the Anti-V5-tag Beads, then screw the top cap on tightly. Tap the tube gently several times. Incubate for 5 minutes at room temperature. Remove the top cap and bottom plug. Centrifuge for 10 seconds.

14. For the second elution, it is not necessary to place the bottom plugs and top cap on the Spin Column. Add 20 μ L Elution Peptide Solution to the Anti-V5-tag Beads, then tap the tube gently several times. Incubate for 1 minute at room temperature. Centrifuge for 10 seconds. The two eluates (step 13 and 14) may be pooled in one sampling tube.


(Note: Steps 11-14 of this protocol are identical to steps 15-18 of the first protocol.)

Related Products:

3317A	V5-tagged Protein Purification Kit Ver.2	2 purifications
3318	V5-tagged Protein Purification Gel Ver.2	1 mL x 1
3315-205	V5-tag peptide	2 mg x 5
3305	c-Myc-tagged Protein Mild Purification Kit ver.2	20 purifications
3305A	c-Myc-tagged Protein Mild Purification Kit ver.2	2 purifications
3306	c-Myc-tagged Protein Mild Purification Gel with Elution Peptide	
	Gel: 1 mL x 1, Peptide: 1 mg x 1	
3307	c-Myc-tagged Protein Mild Purification Gel with Elution Peptide	
	Gel: 1 mL x 5, Peptide: 1 mg x 5	
3300-205	c-Myc-tag peptide	1 mg x 5
3306K	c-Myc-tagged Protein PURIFICATION CARTRIDGE	1 mL x 1
3310	His-tagged Protein Purification Kit	20 purifications
3310A	His-tagged Protein Purification Kit	2 purifications
3311	His-tagged Protein Purification Gel with Elution Peptide	
	Gel: 1 mL x 1, Peptide: 2 mg x 5	
3312	His-tagged Protein Purification Gel with Elution Peptide	
	Gel: 1 mL x 5, Peptide: 2 mg x 25	
3310-205	His-tag peptide	2 mg x 5
3320	HA-tagged Protein Purification Kit	20 purifications
3320A	HA-tagged Protein Purification Kit	2 purifications
3321	HA-tagged Protein Purification Gel	1 mL
3320-205	HA-tag peptide	2 mg x 5
3325	DDDDK-tagged Protein Purification Kit	20 purifications
3325A	DDDDK-tagged Protein Purification Kit	2 purifications
3326	DDDDK-tagged Protein Purification Gel with Elution Peptide	
	Gel: 1 mL x 1, Peptide: 1 mg x 5	
3327	DDDDK-tagged Protein Purification Gel with Elution Peptide	
	Gel: 1 mL x 5, Peptide: 1 mg x 25	
3328R	DDDDK-tagged Protein Purification Gel	5 mL x 1
3329R	DDDDK-tagged Protein Purification Gel	25 mL x 1
3325-205	DDDDK-tag peptide	1 mg x 5
3326K	DDDDK-tagged Protein PURIFICATION CARTRIDGE	1 mL x 1
M215-3	Anti-V5-tag mAb (OZA3)	
M215-7	Anti-V5-tag mAb-HRP-DirecT (OZA3)	

Example of Purification Results 1

Purification of C-terminus V5-tagged TPO protein from culture supernatant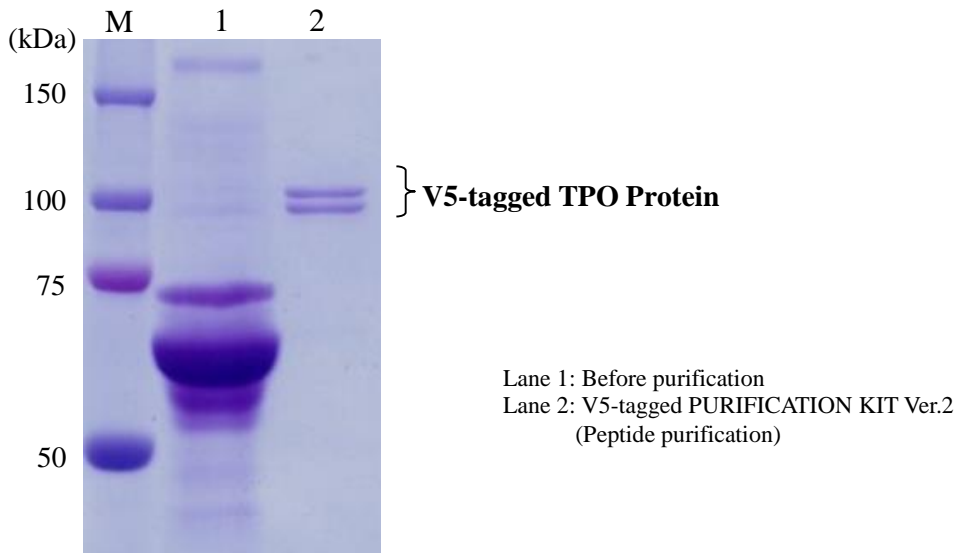


SDS-PAGE (Coomassie Brilliant Blue Staining)

Stable transfectant of *Drosophila melanogaster* Schneider No. 2 cells expressing C-terminus V5-tagged TPO protein were cultured for 3 days in Shields and Sang M3 Insect medium containing 10% Fetal bovine serum. C-terminus V5-tagged TPO protein was purified from 1 mL of cultured medium according to the preceding protocol B. For comparison, elution was carried out not only with peptide solution but also with acid solution and alkaline solution. Each purification was conducted with the same amount of Anti-V5-tag Beads (5 μ L) and the same amount of elution solution (20 μ L x 2 times and then pooled).

Example of Purification Results 2

Purification and enzymatic activity of N-terminus V5-tagged β -galactosidase

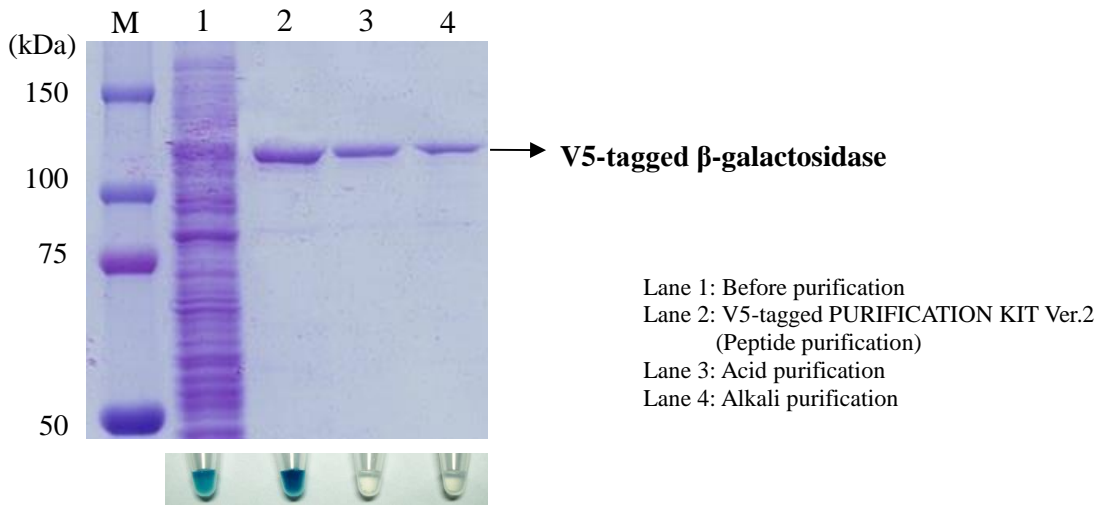SDS-PAGE (Coomassie Brilliant Blue Staining) & X-gal staining

Human embryonic kidney cells (293T) were transfected with pcDNA-V5- β -galactosidase and cultured for 60 hours. Cells were then lysed in the Lysis buffer (1 mL/100-mm dish) and purified according to the preceding protocol A. For comparison, elution was carried out not only with peptide solution but also with acid solution and alkaline solution. Each purification was conducted with the same amount of Anti-V5-tag Beads (5 μ L) and the same amount of elution solution (20 μ L x 2 times and then pooled).

Elution Peptide Solution: neutral pH

0.1 M Glycine-HCl: pH 3.0 (Neutralize the elution immediately with 1 M Tris-HCl, pH 8.0)
(Acid elution solution)

0.1 M NH_3 : pH 11.3 (Neutralize the elution immediately with 1 N acetic acid)
(Alkali elution solution)

Enzymatic activity of each purified proteins was performed using standard X-gal staining method.

Additional Information

Several reagents were examined whether or not they were suitable for use with the V5-tagged Protein PURIFICATION KIT Ver.2. For example, RIPA buffer could be used for preparation of cell lysate. The results are listed below.

Chaotropic agents

Urea	1 M	Yes
Guanidine-HCl	1 M	Yes

Reducing agents

DTT	10 mM	Yes
2-Mercaptoethanol	10 mM	Yes

Surfactants

Nonionic	Tween-20	5%	Yes
	Triton X-100	5%	Yes
	NP-40	1%	Yes
	Digitonin	1%	Yes
	n-Octyl-β-D-galactoside	1%	Yes
Zwitterionic	CHAPS	1%	Yes
	CHAPSO	1%	Yes
Anionic	SDS	0.1%	Yes
	Sodium Deoxycholate	0.5%	Yes

Others

NaCl	1 M	Yes
Glycerol	10%	Yes
EDTA	10 mM	Yes

These reagents can be used in the Lysis buffer for this kit, up to the indicated concentration.

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3317A	V5-tagged Protein Purification Kit Ver.2	2 purifications
3318	V5-tagged Protein Purification Gel Ver.2	1 mL x 1
3315-205	V5-tag peptide	2 mg x 5
3305	c-Myc-tagged Protein Mild Purification Kit Ver.2	20 purifications
3306	c-Myc-tagged Protein Mild Purification Gel with Elution Peptide	
	Gel: 1 mL x 1, Peptide: 1 mg x 1	
3300-205	c-Myc-tag peptide	1 mg x 5
3306K	c-Myc-tagged Protein PURIFICATION CARTRIDGE	1 mL x 1
3310	His-tagged Protein Purification Kit	20 purifications
3311	His-tagged Protein Purification Gel with Elution Peptide	
	Gel: 1 mL x 1, Peptide: 2 mg x 5	
3310-205	His-tag peptide	2 mg x 5
3320	HA-tagged Protein Purification Kit	20 purifications
3321	HA-tagged Protein Purification Gel	1 mL
3320-205	HA-tag peptide	2 mg x 5
3325	DDDDK-tagged Protein Purification Kit	20 purifications
3326	DDDDK-tagged Protein Purification Gel with Elution Peptide	
	Gel: 1 mL x 1, Peptide: 1 mg x 5	
3328R	DDDDK-tagged Protein Purification Gel	5 mL x 1
3329R	DDDDK-tagged Protein Purification Gel	25 mL x 1
3325-205	DDDDK-tag peptide	1 mg x 5
3326K	DDDDK-tagged Protein PURIFICATION CARTRIDGE	1 mL x 1
M215-3	Anti-V5-tag mAb (OZA3)	
M215-7	Anti-V5-tag mAb-HRP-DirecT (OZA3)	

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